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E-Mail: info@orsatec.de

<http://www.eukitt.de>

Instructions EUKITT® mounting medium

EUKITT™ is a mounting medium for microscopy and for in vitro diagnostics. It is used for bonding and sample preservation during manual or automated coverslipping and can be used as special adhesive and sealant with glass, non-plastic and other non-reactive (inert) materials. It spreads evenly and without air bubbles.

Characteristics:

Quick-drying, refractive index close to glass (1.490), chemically neutral, low water-sensitive, no reaction to staining, temperature range: 42 °C - 17 °C, resistant to UV light, yellowing-resistant, cold-resistant, optically clear for 10 years, used in place of epoxy resins during mounting of petrographic thin sections, as with fluorescent dyes in bone. Store at room temperature.

Safety Instructions:

The diagnostic application of EUKITT® and the evaluation of the preparations carried out by a physician or according to suitably qualified person are the responsibility of the laboratory doctor. The respective national legislation must be observed. Likewise, accident prevention and hygiene regulations must be observed.

The diagnostic application of EUKITT® and the evaluation of the preparations carried out by a doctor or for qualified person in the responsibility of the laboratory doctor.

National law must be respected. Additionally, accident prevention and hygiene regulations must be respected.

Warning:

Before use, please read all information carefully.

Biohazard:

Pay attention to the use of appropriate personal protective equipment when handling potentially infectious materials.

Do not use, if package is damaged:

Damage of the packaging can lead to the emission of EUKITT®. Take attention to the risk of wetting and take appropriate safety measures (for example, wearing hand gloves).

Usability:

The usability is limited to 36 months. Expired must not be used any longer for diagnostics.

Instructions:

EUKITT® is ready to use. After adaptation to room temperature it can be used without further preparation steps.

Pay attention to instructions:

The use of EUKITT® should be carried out according to the use instructions of the respective system or reagent supplier or to your own validated procedures.

-Refractive index $n[20\text{ °C}]$: 1.49

-Viscosity: 250-450 mPa*s

-Drying time: 20 minutes

Warning:

Only properly trained persons are permitted to use EUKITT®.

Since mounting media like EUKITT® are not used by amateurs or for self-testing, we dispense with the description of the many ways of usage and refer to the relevant technical and training literature.

Storage instructions:

Optimally, EUKITT® is stored at 20 °C. After use, it is advisable to close the bottle immediately. In regions of high temperatures thickenings can occur due to evaporation of the solvent, which can easily be remedied by dilution with xylene. Required temperature changes should be carried out slowly. Due to the limited period of use the older stock items should always be used first.

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Flammable liquid and vapour (H226). Harmful in
contact with skin (H312). Harmful if inhaled (H332).
Causes skin irritation (H315). Causes serious eye
irritation (H319). May cause respiratory irritation
(H335). May cause damage to organs through
prolonged or repeated exposure (H373).
Keep away from heat/sparks/open flames/hot sur-
faces. No smoking (P210). Avoid breathing vapours
(P261). Wear protective gloves /protective clothing/
eye protection / face protection (P280). IF ON SKIN
(or hair): Remove/Take off immediately all contami-
nated clothing. Rinse skin with water/shower
(P303+P361+P353). Dispose of contents/container
in accordance with local regulation (P501). Contains
methyl methacrylate, n-butyl methacrylate. May
produce an allergic reaction (EUH208).

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Flammable liquid and vapour (H226). Causes skin irritation (H315). May cause an allergic skin reaction (H317). Very toxic to aquatic life with long lasting effects (H410).

Keep away from heat/sparks/open flames/hot surfaces. No smoking (P210). Avoid breathing vapours (P261). Wear protective gloves /protective clothing/ eye protection / face protection (P280). IF ON SKIN (or hair): Remove/Take off immediately all contaminated clothing. Rinse skin with water/shower (P303+P361+P353). Store in a well-ventilated place. Keep cool (P403+P235). Dispose of contents/container in accordance with local regulation (P501).

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